**Signs**  
Included in the packaging of the BROEN-LAB emergency shower is the customer receives an international emergency symbol.  
This symbol must be placed where it is visible and where it clearly indicates the installation of an emergency shower.

**Instructions for Use**

- Start flushing immediately - the second count
- Remove contaminated clothing which is wet and dark to the skin
- Continuous flushing
- In case of an eye injury keep the eyelids open when continuing flushing
- Call for an ambulance. Advise the doctor the accident
- Continuous flushing in the ambulance
- Keep the flushing unit emergency personnel during transport to the special ward

**Do remember to inform all employees about:**

- How to prevent accidents
- How to react in an emergency
- The location of the emergency shower
- How to operate the emergency shower
- How to help injured persons in the best way

**Warning!**  
Before the installation of your BROEN-LAB emergency shower, all staff who is responsible for the safety of the product, and the system was found not to have been followed out, the warranty will be void.

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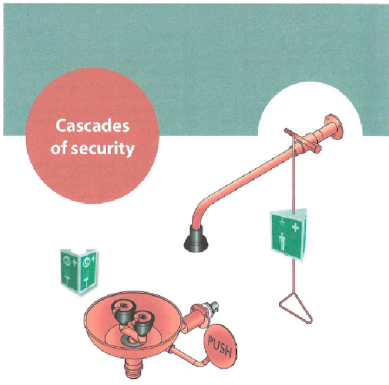
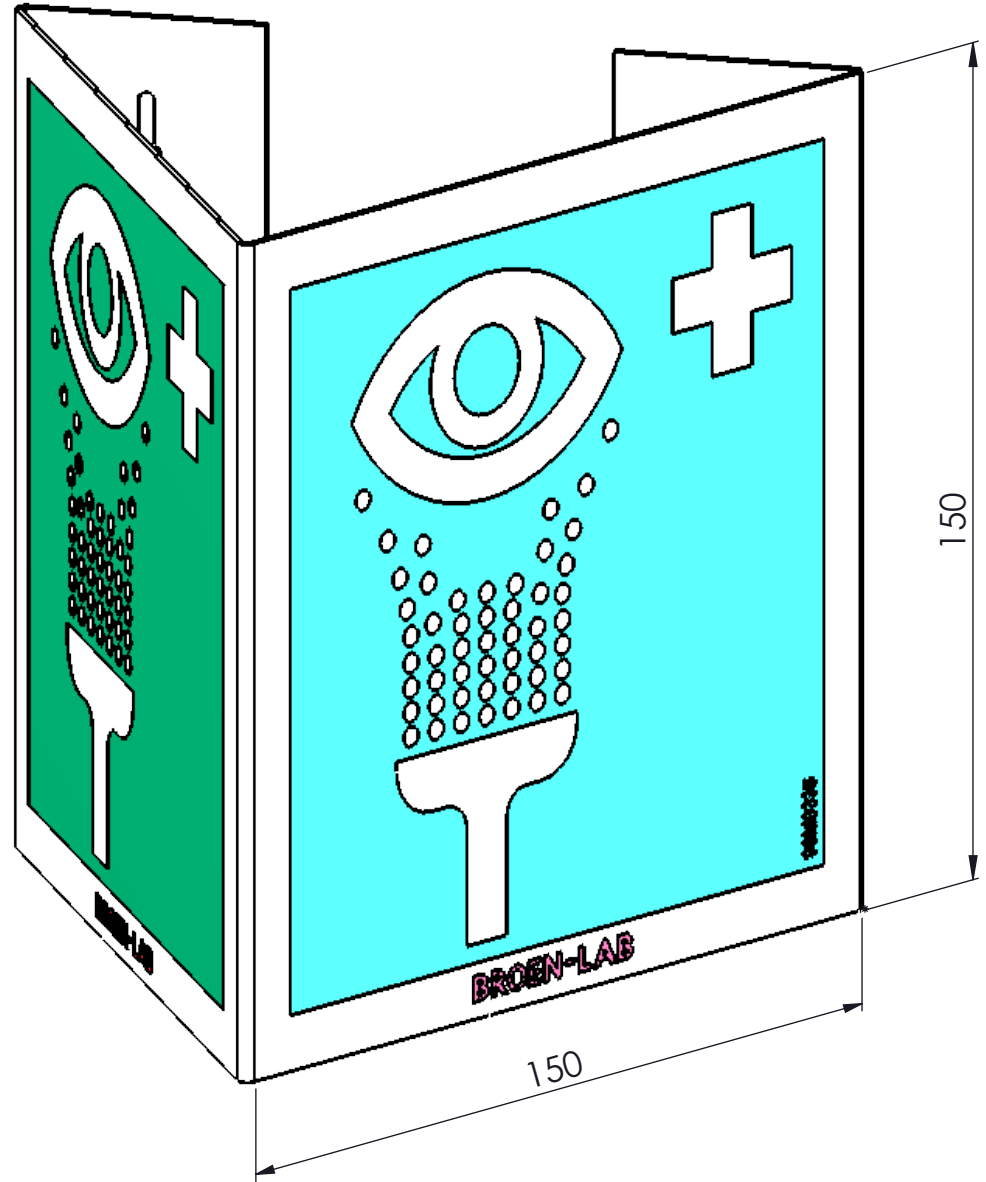
Drawing type: Customer drawing

Weight: Kg

All dimensions marked with  are inspection dimensions

Unspecified linear tolerances:

Unspecified angular tolerances:

Designed by: KEC  
Drawn by: HEM

Released on TM No.

Date: 24-03-20

Format: A3  
Drawing No.:

17001970C

Rev.	Pos.	Date	TM No.	Design by	Drawn by	Change:
A		30-01-07	5086	Admin	Updatet signs	

**BROEN-LAB**

This drawing is our property and cannot be given to a third party, copied or otherwise used without our permission. BROEN-LAB A/S